



PTR/HT Filters

High Temperature PTFE Membrane



PTR/HT filters are validated Polytetrafluoroethylene (PTFE) membrane cartridge filters assembled with high temperature polypropylene components. They are used for sterilizing in process gas applications and tank vents operated at elevated temperatures (up to 221°F). The PTR/HT membrane has a pore size of 0.22 μm .

These hydrophobic PTR/HT filters have high air flow and low pressure drops. Each cartridge module is individually integrity tested using the water intrusion method before it is released from manufacture.

Critical Process provides unrivaled delivery times, technical consulting before purchasing, and very competitively priced high-performance products. Our comprehensive testing & analysis and validation services support your team whenever they need it. Your process experts partnering with our filtration experts is how we deliver your company's solution right the first time.

Sterilizing Filters


Tank Vent & Process Gas



CARTRIDGES – Nominal Dimensions

Length: 5 to 40 in. (12.7 to 101.6 cm)

Outside Diameter: 2.75 in. (7.0 cm)

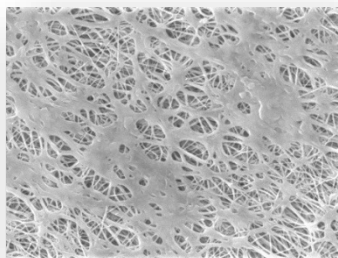

PTR/HT filters are recommended for:

- Tank Vents
- Compressed Air
- Pressurized Gases
- Fermentation Air

Operated at elevated temperatures

Maximum Operating Parameters

CARTRIDGES	
Operating Temperature	221 °F (105 °C)
Forward Differential Pressure	80 psid at 68 °F (5.52 bard at 20 °C)
Reverse Differential Pressure	50 psid at 68 °F (3.45 bard at 20 °C)
Recommended Maximum Service Life	1 Year

Sanitization & Sterilization

Inline Steam	275 °F (135 °C), 30 min, 25+ cycles
Autoclave	250 °F (121 °C), 30 min, 25+ cycles
Chemical Sanitization	Performed using industry standard concentrations of hydrogen peroxide, peracetic acid, sodium hypochlorite and other selected chemicals.

Filtration Area (Nominal)

CARTRIDGES					
Length	5"	10"	20"	30"	40"
	12.7cm	25.4cm	50.8cm	76.2cm	101.6cm
Area	3.8 ft²	8.2 ft²	16.4 ft²	24.6 ft²	32.8 ft²
	0.36m²	0.76m²	1.52m²	2.29m²	3.05m²


Integrity Testing

PORE SIZE	WATER INTRUSION TEST PRESSURE		BUBBLE POINT MINIMUM*	
	PSIG	BARG	PSIG	BARG
0.22	35	2.41	18	1.24

WATER INTRUSION SPECIFICATIONS (mL/10 min)					
Length	5"	10"	20"	30"	40"
0.22µm	≤ 5.6	≤ 13	≤ 26	≤ 39	≤ 52


* Bubble Point for membrane wetted with 60% IPA / 40% water

Water Flow Rates for PTR/HT Cartridges by Pore Size

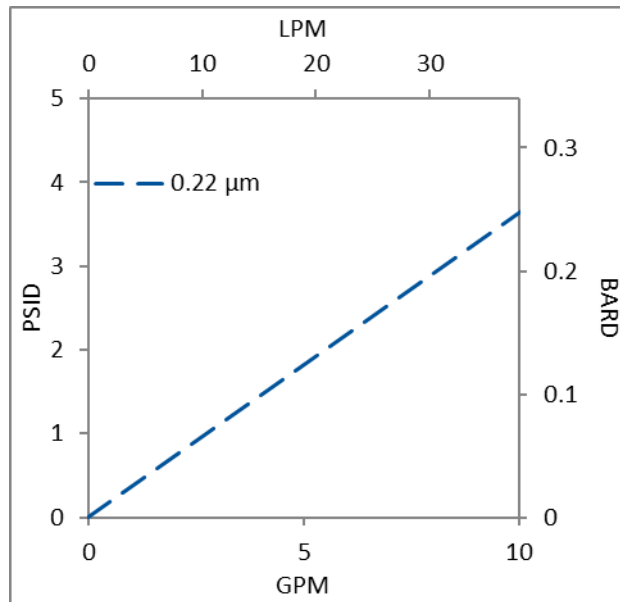

Flow rates for Cartridge filters are per 10-inch length. The test fluid is water at ambient temperature.

Air/Gas Flow Rates for PTR/HT Cartridges by Pore Size

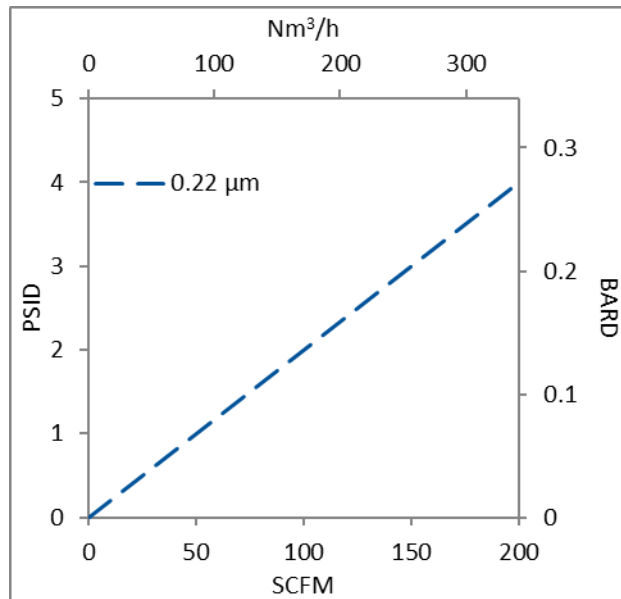

Flow rates for Cartridge filters are per 10-inch length. The test fluid is compressed air at ambient temperature.

Construction Materials

Filtration Media	Polytetrafluoroethylene (PTFE) Membrane
Media Support	High Temperature Polypropylene
End Caps, Center Core, Outer Support Cage	High Temperature Polypropylene
Sealing Method	Thermal Bonding
O-Rings/Gaskets Cartridges only	Buna, Viton® (or FKM), EPDM, Silicone, FEP Encapsulated Silicone, FEP Encapsulated Viton (or FKM)

Validation

PTR/HT cartridges are validated using test procedures that comply with the intent of BFE protocols for the determination of bacterial retention in filters used for air and gas filtration. The challenge level for the 0.22 µm filter membrane is 7.5×10^7 organisms per cm² of filter media of *Brevundimonas diminuta* (ATCC 19146).

PTR/HT filters are also validated using test procedures that comply with ASTM F 838-15(ae1) protocols for the determination of bacterial retention in filters used for liquid filtration. The challenge level is a minimum of 10^7 *Brevundimonas diminuta* organisms per cm² of filter media. CPF filters have > 7-log removal when challenged with the *Brevundimonas diminuta* (0.22µm meets the FDA definition of sterilizing grade filters).

Validation Guides available upon request.

Endotoxins

The levels of bacterial endotoxins in aqueous extracts from PTR/HT filters are below current USP limits as specified for water for injection.

Extractables

PTR/HT filters typically exhibit low levels of non-volatile residues.

TOC and Conductivity

The PTR/HT filters conform with TOC standards of USP <643> and the water conductivity standards of USP <645> after an appropriate flush with purified water.

Toxicity Compliance

Materials used to construct the PTR/HT filters are non-toxic and meet the requirements for the MEM Elution Cytotoxicity Test and the requirements for Biological Reactivity Tests in the current version of the United States Pharmacopeia (USP) for Class VI - 121 °C Plastics.

Non-Fiber Releasing

PTR/HT filters comply with Title 21 CFR sections 210.3 (b)(6) and 211.72, for non-fiber releasing filters.

FDA Compliance

Materials meet the requirements listed by the FDA as appropriate for use in articles intended for repeated food contact as specified in Title 21 CFR sections 174.5, 177.1500, 177.1520, 177.1630, 177.2440, and 177.2600 as applicable.

PTR/HT Filters Ordering Information

Fill in the corresponding codes in the boxes below to build your Part Number.

To consult with one of our technical team members, request a quote or place an order:
call (603) 880-4420 or [contact us here](#).

Cartridge Filters

PTR		S	OHT			
Pore Size Code -20 = 0.22 μm	Length 05 = 5 in. (12.7 cm) 01 = 10 in. (25.4 cm) 02 = 20 in. (50.8 cm) 03 = 30 in. (76.2 cm) 04 = 40 in. (101.6 cm)	O-Ring/Gasket Code S = Silicone B = Buna V = Viton (or FKM) T = FEP Encapsulated Viton (or FKM) E = EPDM R = FEP Encapsulated Silicone		End Cap Code 0 = Flat Gasket, DOE 2 = 2-222 O-ring/Plug 5 = 2-222 O-ring/Flat 6 = 2-226 O-ring/Flat 8 = 2-222 O-ring/Spear 9 = 2-226 O-ring/Spear		



One Chestnut Street
Nashua, NH 03060
603.880.4420
FAX: 603.880.4536

CriticalProcess.com

The information contained herein is subject to change without notice. The Critical Process Filtration logo is a trademark of Critical Process Filtration, Inc. Viton is a trademark of DuPont Performance Elastomers L.L.C.
© 2020 Critical Process Filtration, Inc. • All Rights Reserved

Data Sheet PTRHTDS Rev A